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LuciSwab™ 3.2-400 Instruction Manual

Product code:60344

CAUTION !
Please make sure to read the precautions and instructions in this Instruction Manual before attempting to use the product and exercise extreme caution when using it.

"LuciSwab 3.2-400" is the long stem cotton swab which is exclusive use for the hygiene monitoring using LuciPac A3 Surface, LuciPac A3 Water, LuciPac Pen and LuciPac Pen-AQUA.

The size of the long stem swab: Cotton bud 3.2mmΦ and stem length 400 mm

[Applications]

This product is suitable for swabbing inside of the channels of medical endoscopes and of filling nozzles in beverage industry, which have not accessible by LuciPac and/or other ATP swab test kits.

[Contents]

The LuciSwab 3.2-400 contains 5 bags of 20 individually packed long stem swabs in one bag.(for a total of 100 swabs)

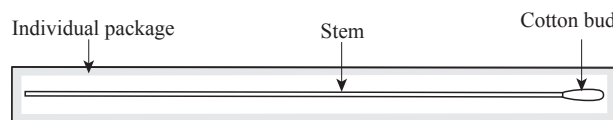
Fig. The name of each part of the long stem swab.

[Precautions for Use]

Please make sure to follow the precautions listed below in order to obtain optimal performance from this product.

- ① LuciSwab may be kept at room temperature.
- ② LuciSwab is specially manufactured for cleanliness testing. Other commercial cotton swabs may not produce correct results.
- ③ Choose LuciSwab 2.8-400 or LuciSwab 3.2-400, depending on the inner diameter of the object to be tested.
- ④ Do not push the LuciSwab into the narrow inspection point forcibly. Stop pushing it forcibly when there is strong resistance at the time of the insertion. If not done, the cotton bud may fall off when you pull the swab up.
- ⑤ Be sure to use only designated products when taking measurements for amounts of luminescence. This product cannot be used with nondesignated products.

[Designated Product and kit for Measurement]

Designated Product : Lumitester PD-20 and Lumitester PD-30

Designated kit: LuciPac A3 Surface , LuciPac A3 Water,

LuciPac Pen , LuciPac Pen-AQUA

(Manufacturer: Kikkoman Biochemifa Company)

*Be sure to use only designated products and kit when taking measurements.

[Disposal Methods]

This product can be disposed of as regular garbage, but when disposing of it, it would be better to separate the parts and dispose of each one properly in accordance with the local regulations outlined by the local governments for proper disposal of waste materials. Moreover, when the product was used to test medical devices, dispose it as clinical waste if needed.

The main materials and parts used in this product are listed below. No PVC materials are used in the making of the plastics in this kit.

Table. Main raw materials of the structural parts of this product

	Structural parts	Raw materials
Long stem swab	Stem	Polyethylene
	Cotton bud	Cotton
	Individual package	Paper, Polyester, Polypropylene
	Outer package	Polyethylene
External packaging	Outer case	Paper
	Shrink-wrap	Polypropylene

[Precautions for Handling]

- ① Exercise caution when storing and disposing of the swab to ensure that none of the substances mix with food or other products.
- ② Please make sure to store this product and its parts out of the reach of young children.

[Warranty]

Kikkoman Biochemifa Company warrants the products in this kit to have a certain level of quality. This warranty guarantees that Kikkoman Biochemifa Company shall replace defective products should any be found. This warranty does not provide any other guarantees. Kikkoman Biochemifa Company shall not be liable for any damages, including special or consequential damages, or expenses arising directly or indirectly from the use of this product.

Symbols used in the packaging and labeling of this product	
	Symbol for "Lot Number." This symbol shall be adjacent to the manufacturer's lot number (e.g. 20160410A) or description of its printed location.
	Symbol for "Consult Instructions Manual."
	Symbol for "Manufacturer." This symbol shall be adjacent to the name and address of the manufacturer.

"LuciPac" and "Lumitester" are registered trademarks of Kikkoman Corporation in Japan and other countries.

Manufacturer

Kikkoman Biochemifa Company

2-1-1 Nishi-shinbashi, Minato-ku, Tokyo 105-0003 , Japan


Phone: +81-3-5521-5492 / Fax: +81-3-5521-5498

E-mail: biochemifa@mail.kikkoman.co.jp


URL: http://biochemifa.kikkoman.co.jp/e/

[Measurement Methods] Note: Please use all LuciPac devices in the same way.


Leave "LuciPac" at room temperature until they have reached room temperature.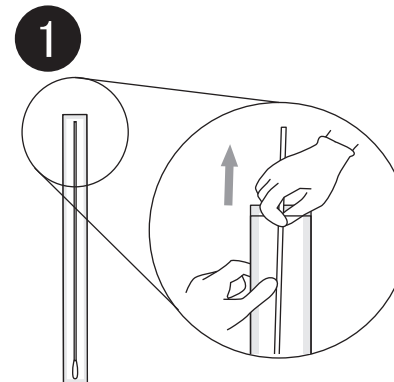


Wear lab gloves to prevent contamination and take out the long stem swab carefully so as to not to touch the parts as this may affect the cleanliness level.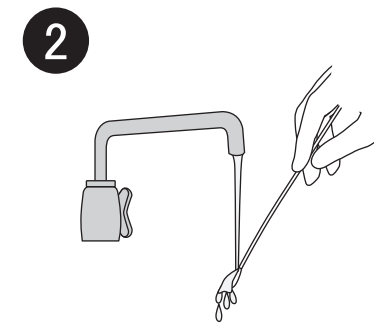


Moisten the cotton bud with tap water. Please avoid using saline solution instead of water.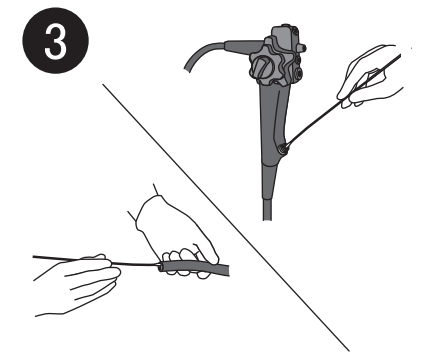


Handle the long stem swab with care so as to not to touch the parts as this may affect the cleanliness level. Insert the long stem swab in the canalicular object to be tested, and wipe it off as long as the length of the long stem swab reaches, and pull the swab up slowly.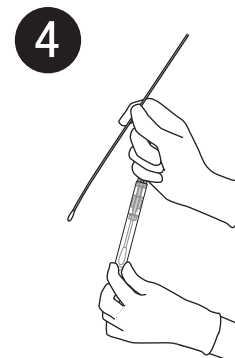


Hold the long stem swab at 12 or 13 cm from the cotton bud, and remove the sampling stick from the main body (casing) of the LuciPac with your free fingers. Handle the sampling stick with care so as not to touch the parts, as this may affect the cleanliness level.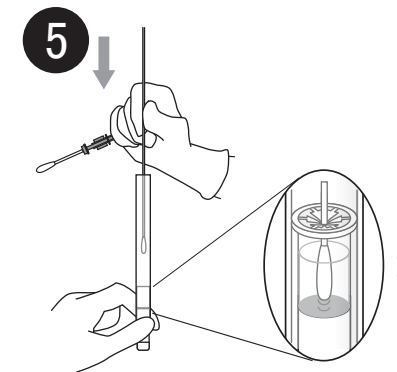

Insert the long stem swab into the main body (casing) of the LuciPac to push and break the center of the upper seal of the container. Be careful not to break bottom seal of the container for releasing reagent. Remove the long stem swab after thoroughly washing the cotton bud in the releasing reagent.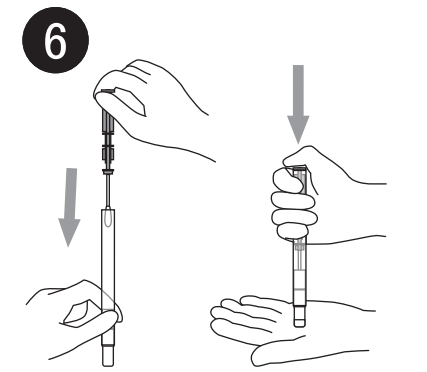

Return the sampling stick into the main body (casing) and push it through all the way by putting the tip of the test tube on the palm of a hand or on a table. (Be careful not to get fingers caught when pushing it.)

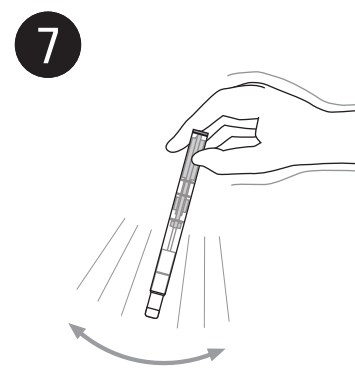Hold the LuciPac casing firmly and shake it. Allow leftover luminescent reagent to thoroughly dissolve.

Insert the LuciPac into the Lumitester. Then, close the Lumitester measurement chamber cover.

Press the ENTER key. Measurement result is displayed after countdown. *Please refer to the instruction manual of LuciPac for data handling.