


## 【USB Driver Installation Procedure】 for Windows10


**Two types of USB drivers** are required to be installed before connecting Lumitester PD-30 and PC.  
Please follow the procedure below.


1 Connect Lumitester PD-30 to PC with the USB cable provided.


### 2 Install the first USB driver.

Click "Start" button and open the "Connected devices".  
Then Click "Device manager".

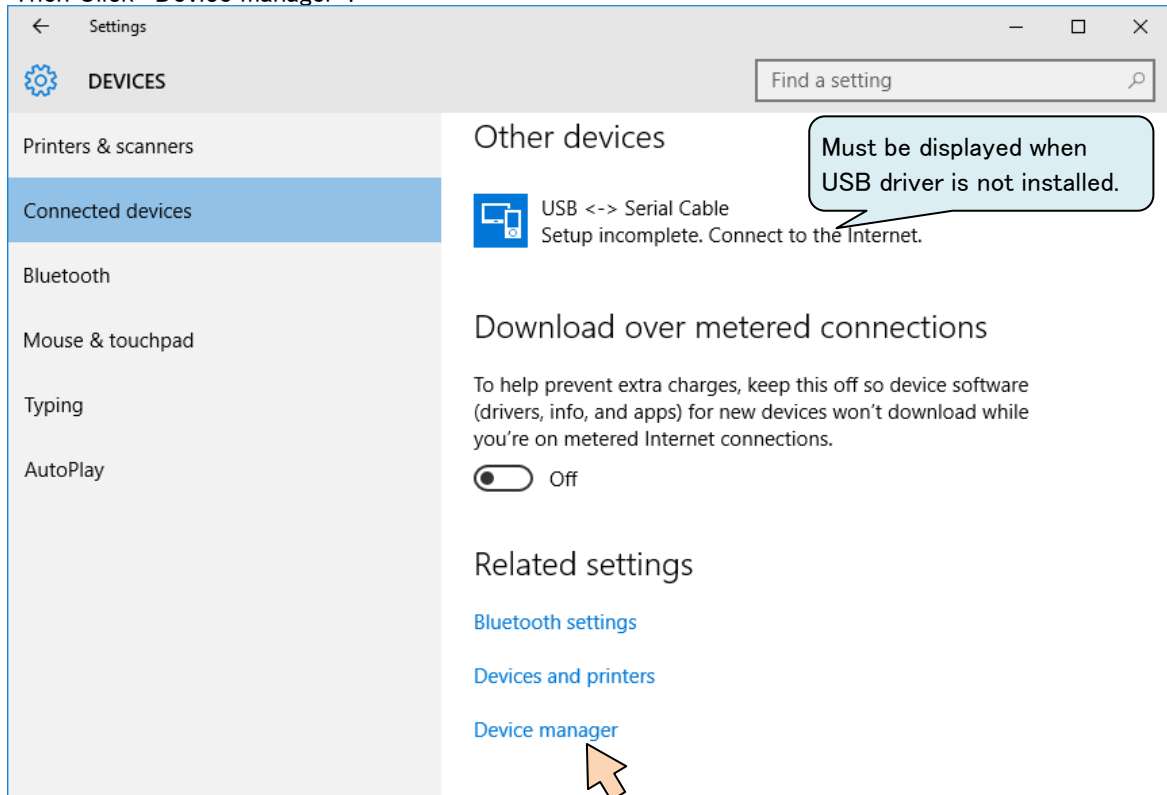


3 The device with the  " will be displayed on Device Manager.


Place the cursor on the name of device with  " and right-click, and select "Update Driver Software".

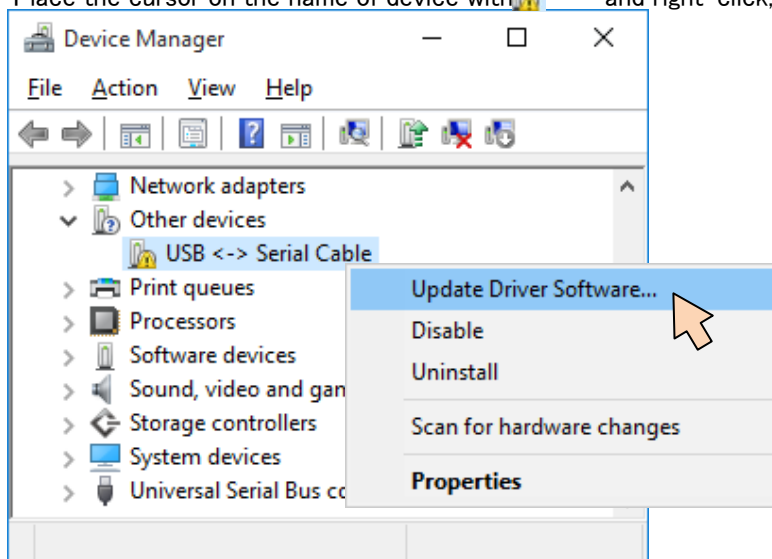


- 4 "Update Driver Software" is displayed on the screen.  
Click "Browse my computer for driver software".

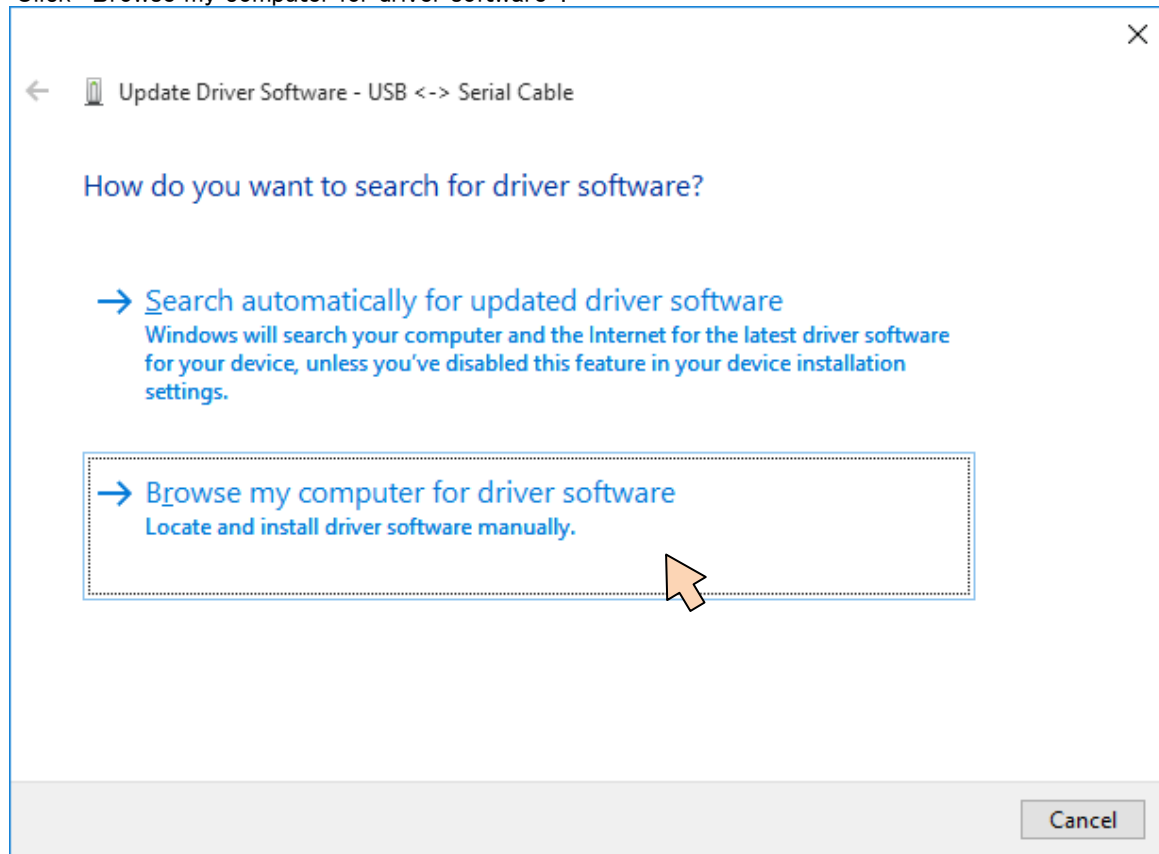

- 5 Click "Browse", find and select the folder of "USB Drivers" downloaded, and then click "Next".

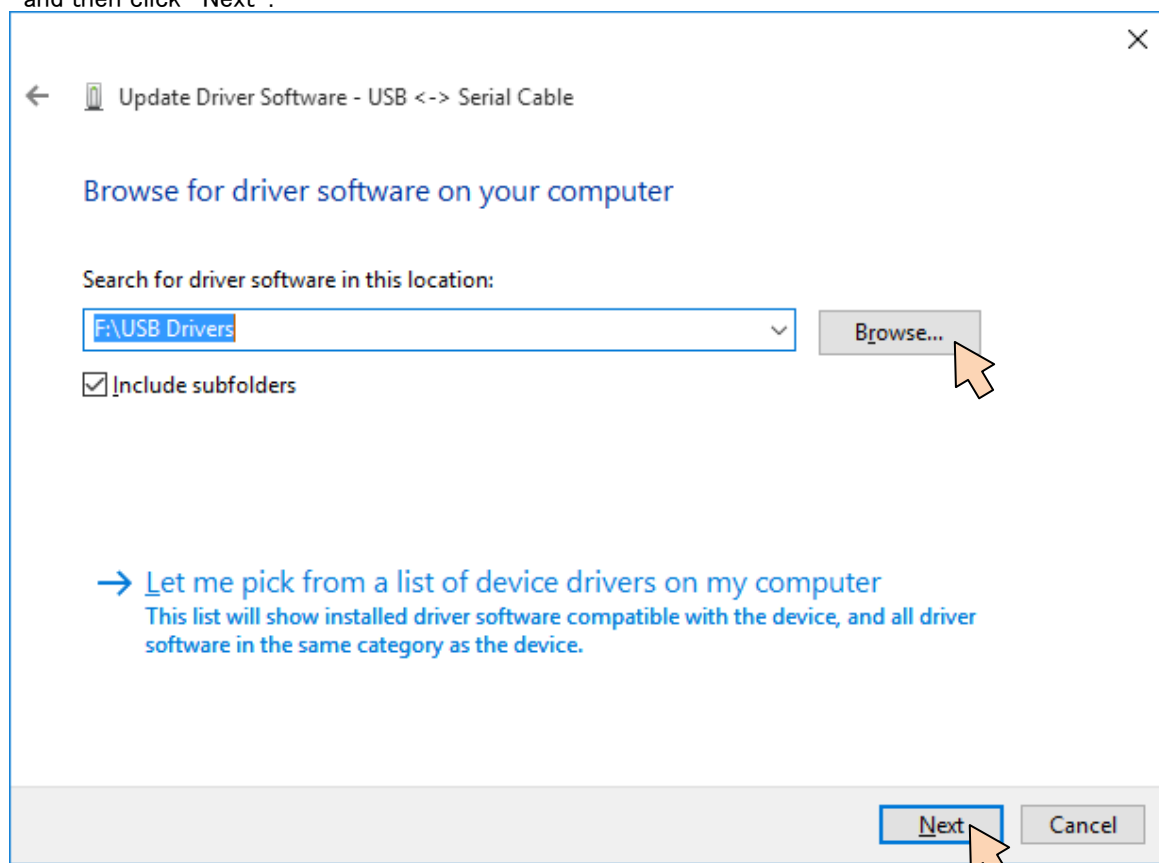


6 Installation of the driver software will begin.

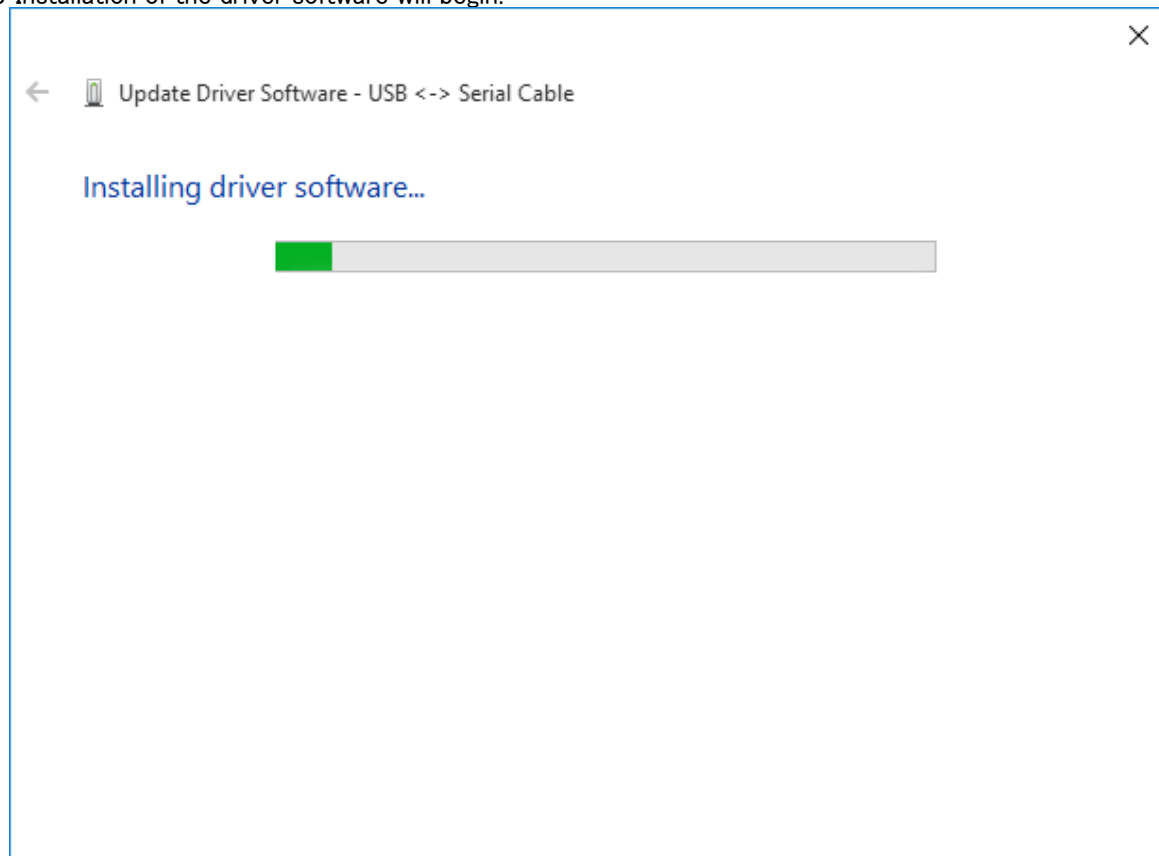


7 The **first USB driver** is now installed.

Click "Close" to close the window. And begin installation for the **second USB driver**.

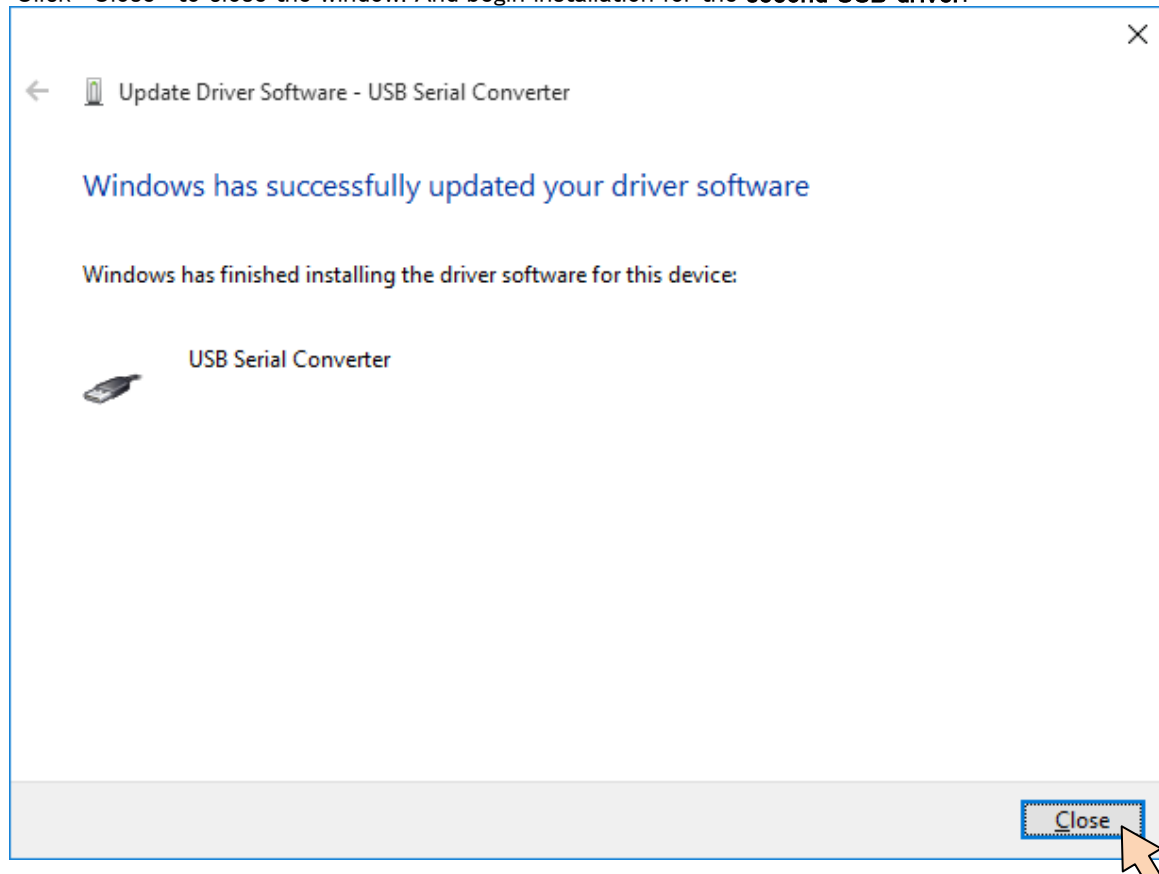

8 For the Installation of **second USB driver**.


Open the Device Manager in the procedure. The device name with the yellow icon is displayed on Device Manager. Place the cursor on the device name and click "Update Driver Software".

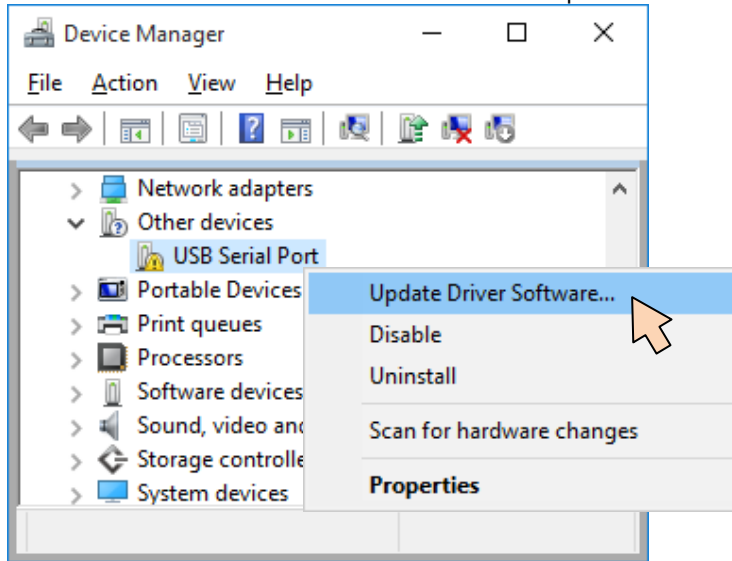

9 Follow the same way as in the procedure 4 to 7 of first USB driver for installation.


When the installation is completed the following screen appears.

Click "Close" to close the window.

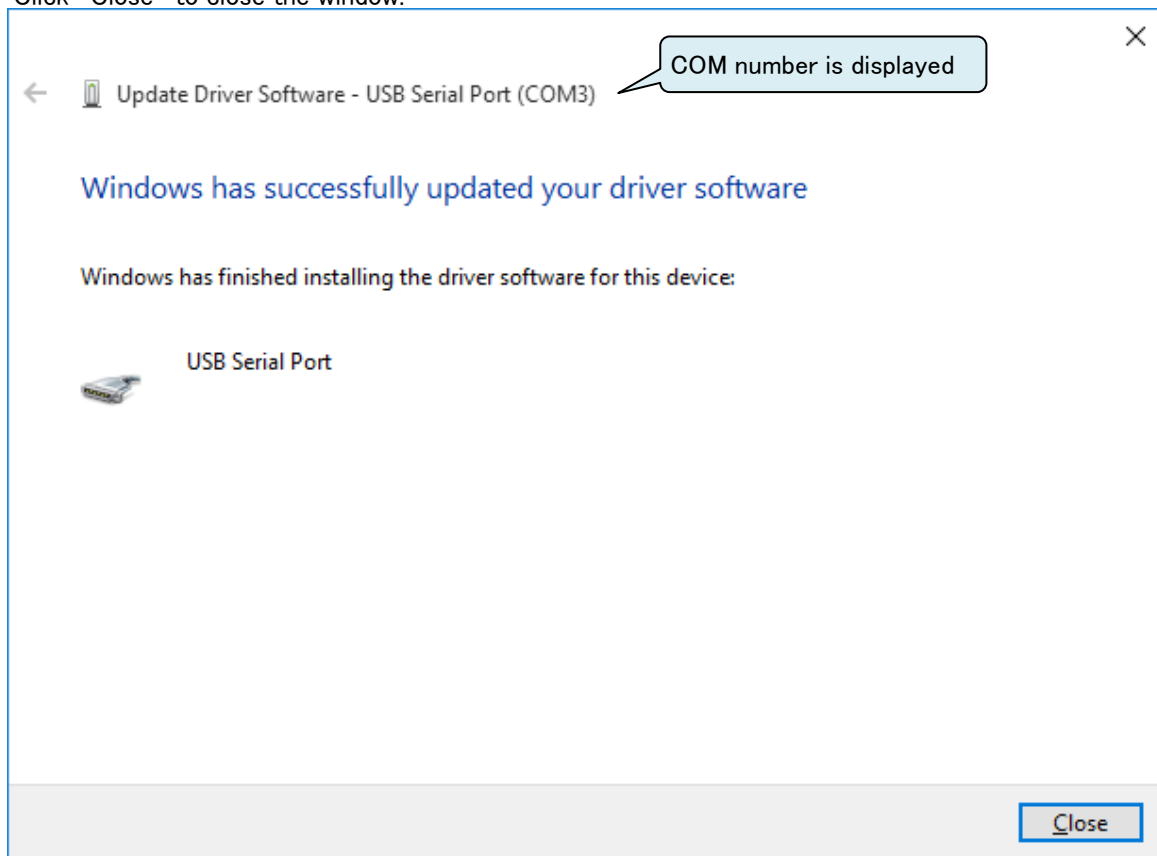


## 【Control Software Installation Procedure】 for Windows10


- 1 Double click "LumitesterPD-30\_Setup.exe" and start.  
Please wait till all the files for Installation is completed.

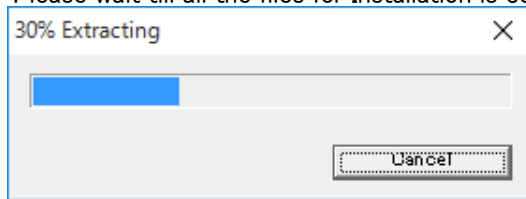


- 2 Click "Lumitester Control Software"

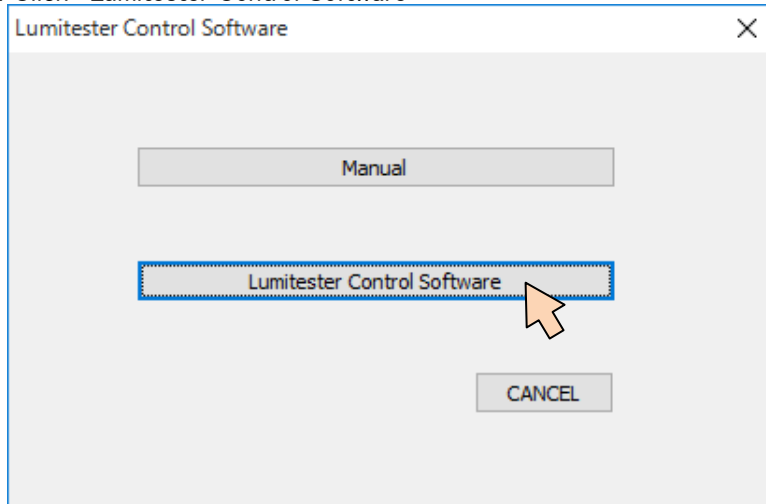


- 3 Click "Yes" when setup is started.

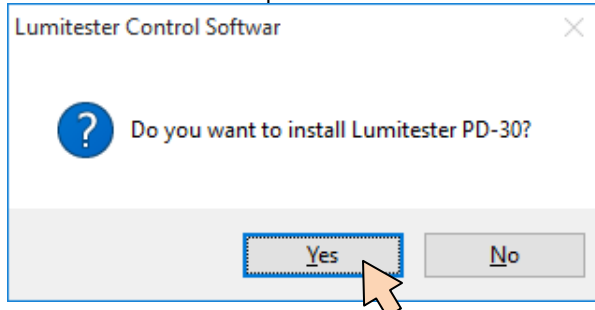


4 Installation is now ready, Click "Next"

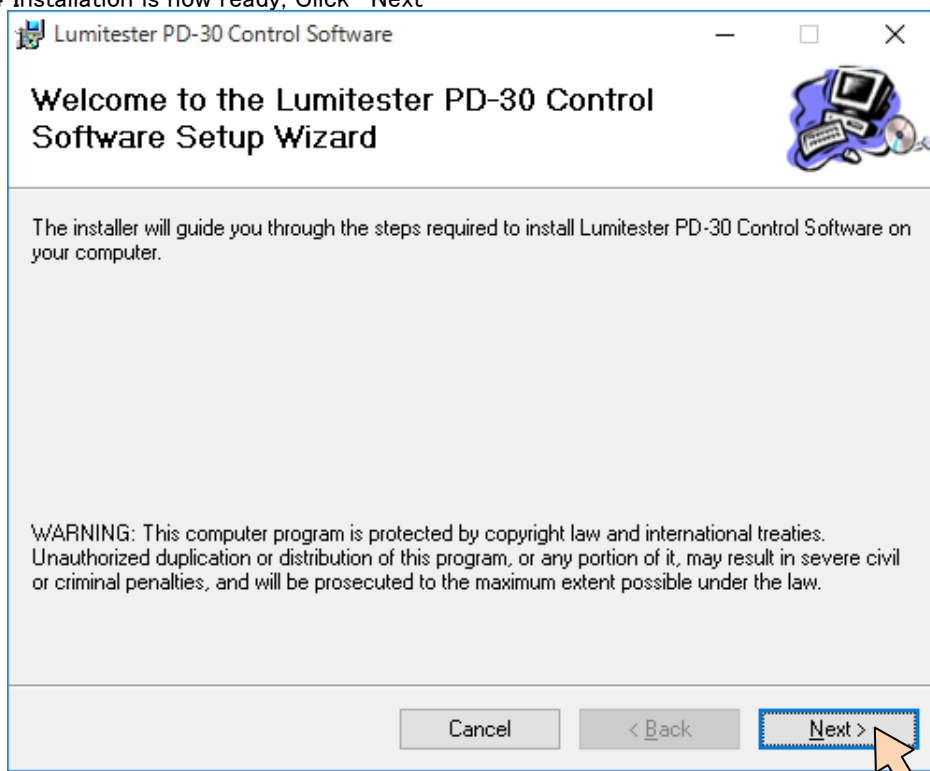


5 Confirm the contents of License.  
Select "Agree" to agree and then click "Next"  
If you do not agree, click "Cancel"

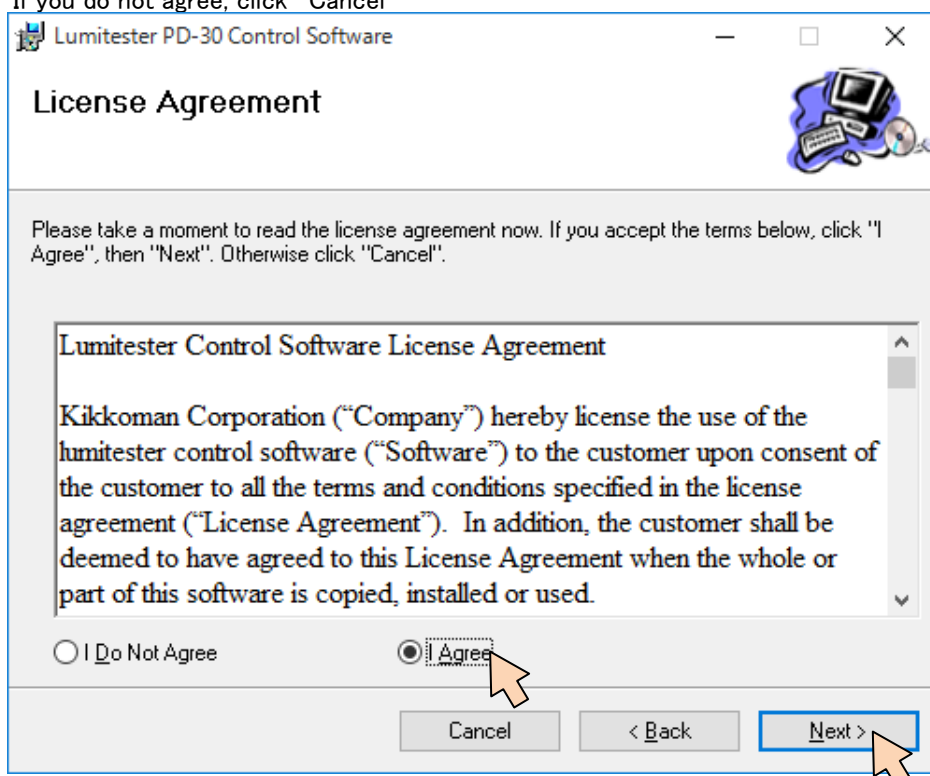


- 6 Click "Yes" if you prefer to have shortcut on the desktop,  
If not, select "No" then click "Next".

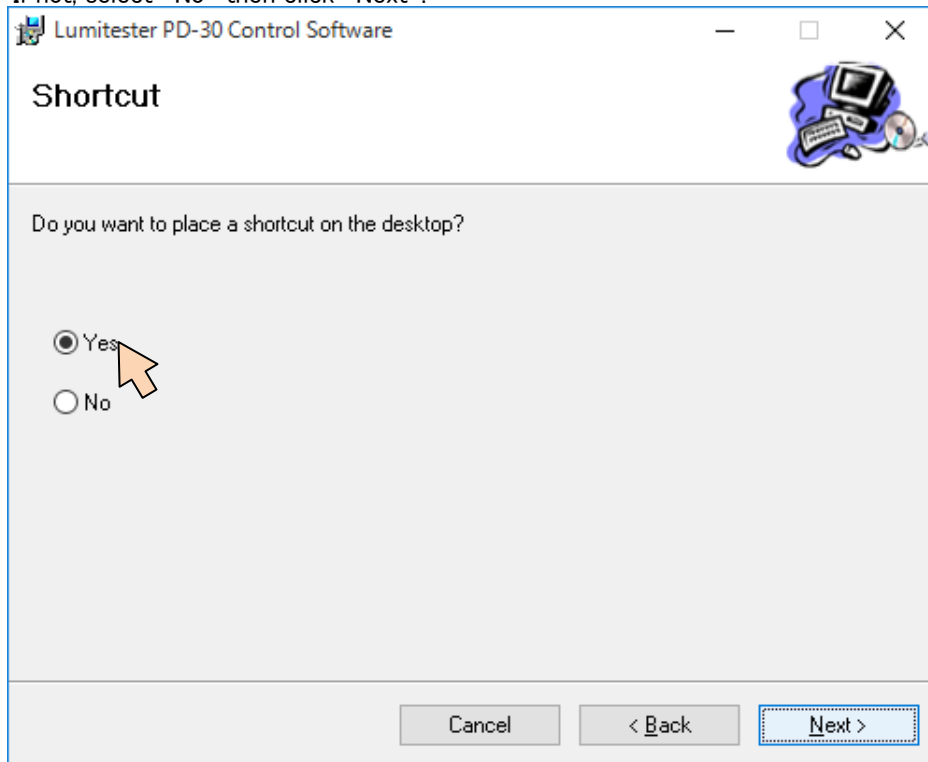


- 7 Enter the name of folder to Install.  
Click "Next" if there are no changes.

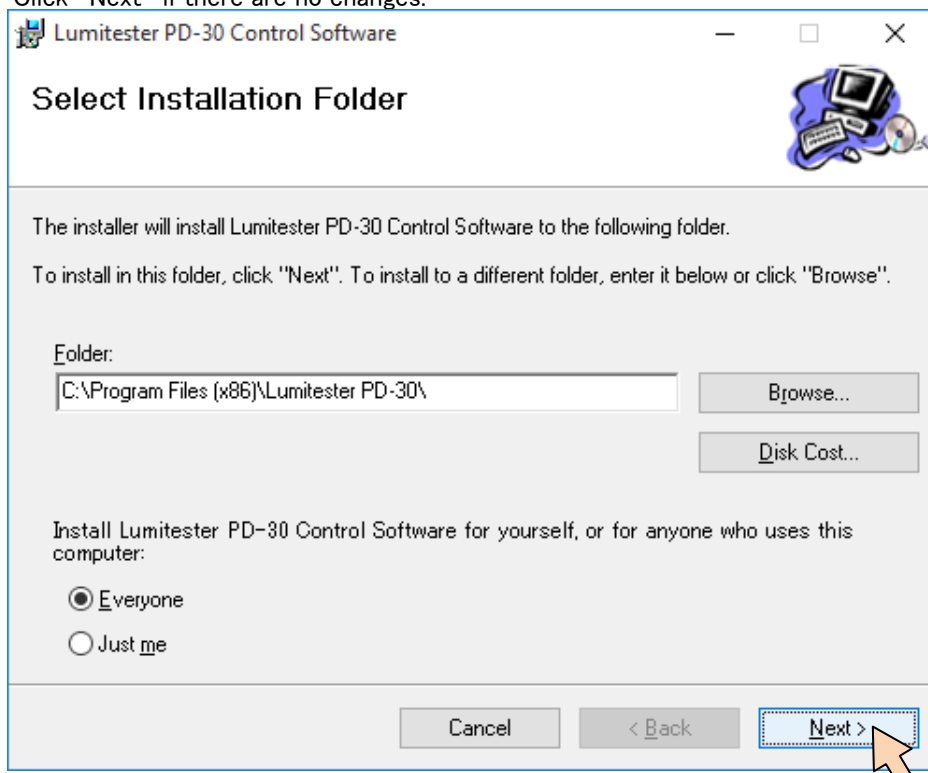

8 Click "Next"

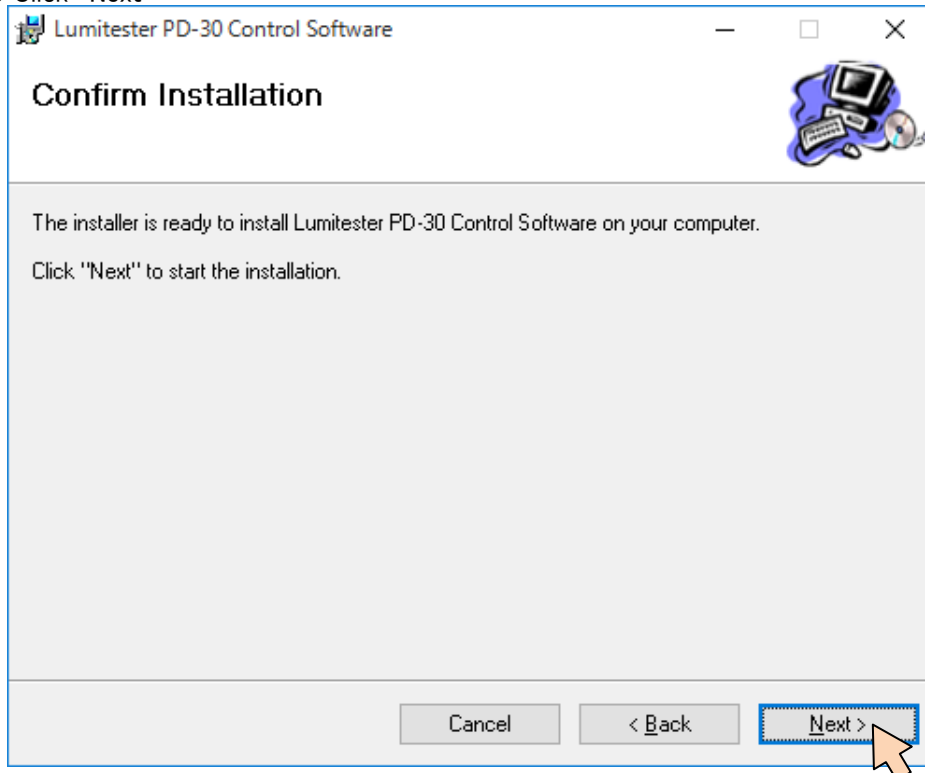

9 Installation is now in progress.

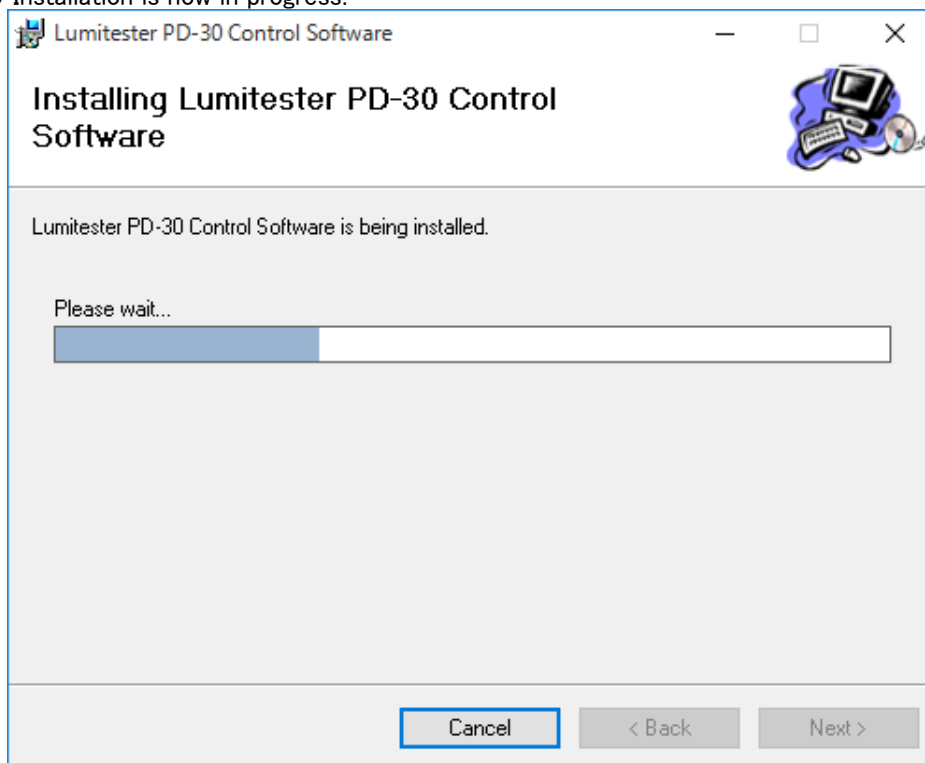



- 10 Installation is complete  
Click "Close" to close the window.

